



# OPERATING INSTRUCTIONS SCREW CAP COMBO VALVE



- TO OPEN VALVE: Use quick opening Screw Cap. Turn cap counter-clockwise less than 1/2 turn (Figure 1). Note: Opening cap more than 1/2 turn may result in leakage.
- TO CONNECT TUBING: Push tubing over end of stem (Figure 2).
- TO CLOSE VALVE: Turn cap clockwise until it stops (Figure 3).
- TO TAKE A SYRINGE SAMPLE: Pierce septum and insert needle approx 3/4" into valve (Figure 4)
- TO REPLACE SEPTUM: Remove screw cap from valve body. Insert a thin, unsharpened rod (I.E. a paper clip) into passageway and push out the orange-colored septum (Figure 5). Insert new septum into the valve cap with the shiny side facing the bag (Figure 6).



FIG. 1 VALVE OPEN



FIG. 2 TUBING CONNECTED



FIG. 3 VALVE CLOSED



FIG. 4 SYRINGE NEEDLE  
IN PASSAGEWAY



FIG. 5 REMOVING  
SEPTUM



FIG. 6 INSERTING  
NEW SEPTUM



1287 Matheson Blvd. E. Tel: 905.629.4545  
Mississauga, Ontario Fax: 905.629.7249  
L4W 1R1 Toll Free: 1.877.427.7837

E-Mail: [info@delta-sci.com](mailto:info@delta-sci.com)  
[www.delta-sci.com](http://www.delta-sci.com)